


# GC CAPILLARY COLUMN APPLICATIONS: ULTRA ALLOY™ STAINLESS STEEL

## Grob Test Mix

COLUMN: UAC-CW, Carbowax  
15M. x 0.25mm I.D. x 0.25µm film  
Cat. No.: UAC-CW-15-0.25F  
Temperature: 60° (3 min. hold) (5°/min.) - 200°C  
Injector: 150°C (on-column)  
Detector: 300°C, FID  
Carrier Gas: 20 ml/min., Helium, U=66cm/sec.

1. dodecane
2. butanediol
3. octanol
4. methyl decanoate
5. dicyclohexylamine
6. methylundecanoate
7. methyldodecanoate
8. 2,6-dimethylaniline
9. 2,6-dimethylphenol
10. ethylhexanoic acid

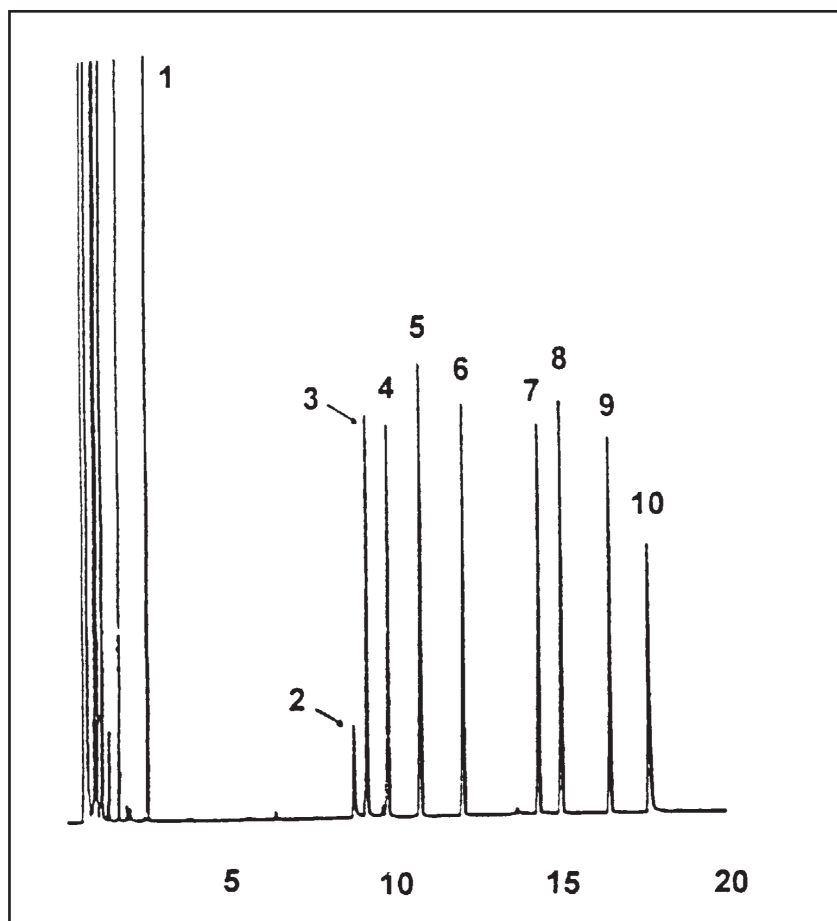

PO Box 3881, Woodbridge, CT 06525 • (203) 393-3112 • FAX (203) 393-0391  
TOLLFREE: 800-275-7033 • E-MAIL: [tech@quadrexcorp.com](mailto:tech@quadrexcorp.com)